## **Declaration of Conformity**

EC-Konformitätserklärung  
EC-Déclaration de conformité  
EC-Dichiarazione di conformità

EC-Declarção de conformidade  
EC-Declaration of conformity  
EC-Declaración de conformidad

**BEL Engineering s.r.l., Via Carlo Carrà 5 Monza 20900 (MB) ITALY**

|           |                                    |                                                                                                                                                     |
|-----------|------------------------------------|-----------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>GB</b> | <b>Declaration of conformity</b>   | <b>We hereby declare that the product to which this declaration refers complies with the following standards.</b>                                   |
| <b>D</b>  | <b>Konformitätserklärung</b>       | <b>Wir erklären hiermit, dass das Produkt, auf das sich diese Erklärung bezieht, mit den nachstehenden Normen übereinstimmt.</b>                    |
| <b>F</b>  | <b>Déclaration de conformité</b>   | <b>Nous déclarons avec cela responsabilité que le produit, auquel se rapporte la présente déclaration, est conforme aux normes citées ci-après.</b> |
| <b>I</b>  | <b>Dichiarazione di conformità</b> | <b>Dichiariamo con ciò che il prodotto al quale la presente dichiarazione si riferisce è conforme alle norme di seguito citate.</b>                 |
| <b>E</b>  | <b>Declarção de conformidade</b>   | <b>Manifestamos en la presente que el product al que se refiere esta declaración està de acuerdo con las normas siguientes..</b>                    |
| <b>P</b>  | <b>Declaración de conformidad</b>  | <b>Declaramos por meio da presente que o produto no qual se refere esta declaração, corresponde às normas seguintes</b>                             |

**Electronic Balance: Series i-Thermo, I-Thermo Touch**

| <b>European Directive</b> | <b>Harmonized Standard</b>             |
|---------------------------|----------------------------------------|
| 2014/30/EU (EMC)          | EN 61326-1:2013                        |
| 2014/35/EU (LVD)          | EN 61010-1:2010<br>EN 61010-2-010:2014 |
| 2011/65/EU (RoHS)         | EN 50581:2012                          |

# QUICK GUIDE

## ELECTRONIC PRECISION MOISTURE BALANCES

Series: i-Thermo L


### ATTENTION:

This is a guide to start quickly working and do NOT substitute the complete user manual; we please you to read completely the user manual for a correct installation and a right use of your new electronic balance.

Usage of the instrument different from the one reported in the user manual does not ensure anymore the safety of the product.

## 1. INSTALLATION INSTRUCTIONS

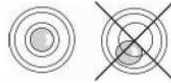- **Remove** the balance, the calibration weight (if included) and the accessories from the package and check if any visible damage of the instrument
- **Do not install** the balance in places with air flows, heavy thermal changes and vibrations. **Place** the support pan and weighing pan on the balance
- **Level the balances** using the levelling feet and looking at the level bubble 
- **Wait for 30 minutes from switch on** and then calibrate the balance (using the appropriate mass if included) following the instructions of the paragraph "calibration"

**WARNING:**

Please follow carefully these steps for installing and use the new balance before starting your work routine. A way of use of the instrument different from this user manual will not guarantee the instrument's safety anymore. Keep this manual in a safe place.

**Please follow these indications to avoid problems and to grant a safe use of moisture analyzer:**

- Use the moisture analyzer exclusively for the moisture determination of samples. Every incorrect use of this instrument may create danger for safety of persons and may cause damages at the instrument or at other objects.
- Do not use the instrument in areas where there is danger of explosions; also make the instrument work only respecting the environment conditions as reported in this manual.
- If this material is used in plants or in environment conditions that require strict safety conditions, please respect the instructions of the directives regarding the installation of this material currently existing in your country.
- This instrument must be used only by qualified personell, who knows the features and characteristics of the sample under test.
- Before starting to work with the instrument for the first time, please check that voltage is the same of your electric mains.
- To get the instrument free from tension, remove the power supply cable.
- Place straight the power supply cable so that to avoid contact with the very hot surfaces of the instrument.
- Use only extension cables that are conform to normatives and that are equipped of a protection conductor.

**Attention, protection against heat**

- Respect the distance and the free space around the instrument to avoid an accumulation of heat in the instrument and the over heating of the instrument itself:
  - 20 cm around the instrument
  - 1 m upon the instrument
- Do not place flammable materials upon, below or near the instrument since the heater warms the surrounding area.
- Remove the samples carefully, the heater and the dishes for samples may be still very hot.

**Danger for persons or objects when working with particular samples: Fire Explosion**

- Flammable or explosive substances
- Substances that contain solvents
- Substances that during drying emit gas or flammable explosive vapors

## 2. Preparing the balance to work

Stainless pan protection  
with ventilation stopper

Underpan

Pan extractor

Pan for the sample

1. Insert stainless pan protection.
2. Insert underpan on the cone of weighth.
3. Insert pan extractor
4. Put the pan for the sample.

1. 15 poles M/F  
cable for  
connection  
balance-heater


2. Balance power supply.  
**Cable VDE**

1. Connect the heater to the balance via cable 15-pin M / F, insert the cable into the two connectors on the rear of the instrument as shown in the figure.
2. Insert the VDE cable in the power connector on the back of the instrument

**N.B. Verify that the power supply indicated on the instrument nameplate matches the one in use in the country where you are installing.**

3. Then connect the VDE cord to the power outlet located near the instrument. Do not use non-compliant with current regulations cables / extensions.


### 3. Keyboard and display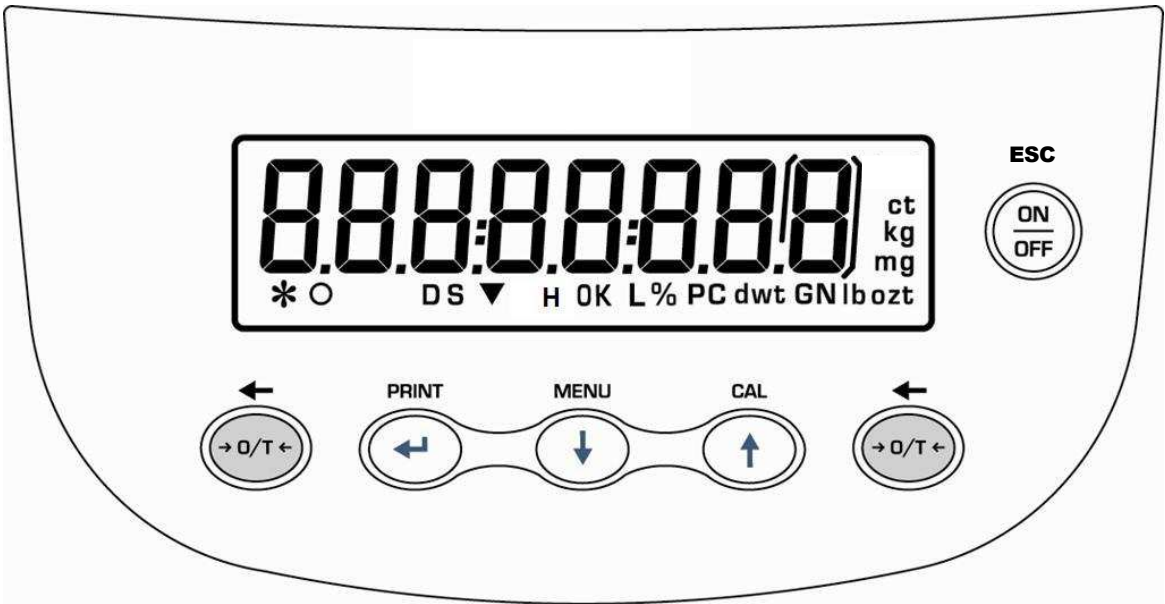


|  |                                                                 |  |                     |
|--|-----------------------------------------------------------------|--|---------------------|
|  | Standby (OFF) or power on (ON) button<br>Escape function button |  | Stability indicator |
|  | TARE or zero button .                                           |  | Zero indicator      |
|  | Selection CONFIRM or SEND data to printer button.               |  | Percentage weighing |
|  | Balance setup MENU button, to set balance's parameters.         |  | Insert data mode    |
|  | Balance CALIBRATION button.                                     |  | Heater is working   |
|  |                                                                 |  | Measure unit        |

## 4. Inputs and outputs

### Rear side and bottom side of the balance model i-Thermo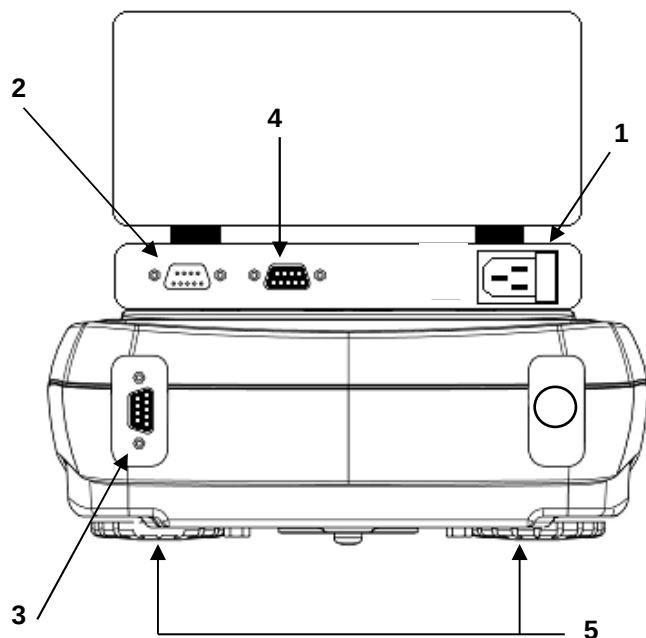


1. **CONNECTOR 1**  
Balance's power supply (VDE)
2. **CONNECTOR 2**  
Connect heater to balance.
3. **CONNECTOR 3**  
Connect balance to heater.
4. **CONNECTOR 4**  
9 pin RS232 Interface for PC and printer
5. **ADJUSTABLE FEET**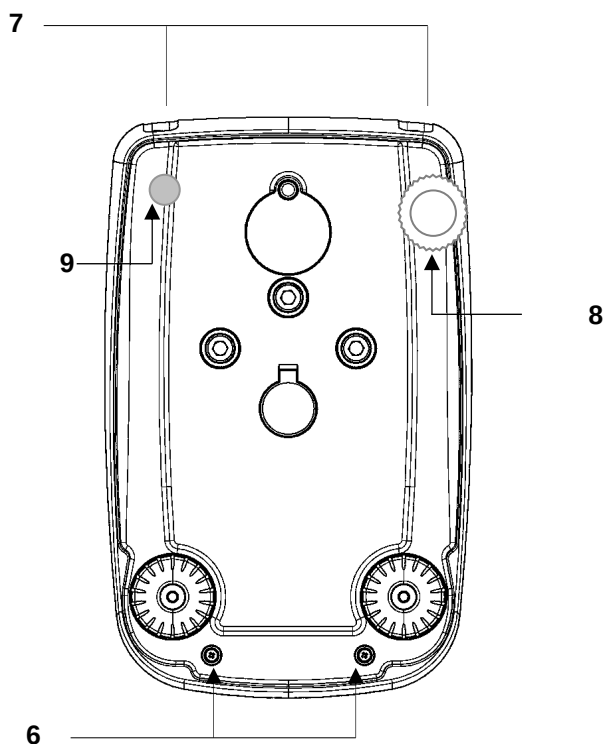

6. **N°2 closing screws of the balance**
7. **N°2 closing screws of the balance: first remove the screws remove adjustabel rear feet(8) and fixed rear feet(9).**
8. **Adjustable rear feet**
9. **Fixed rear feet**

## 5. WEIGHING


In “**STAND BY**” mode press the **ON/OFF** button. Place goods to be weighed on the pan, wait the stability symbol ✱ (asterisk) to appear and then read the weight value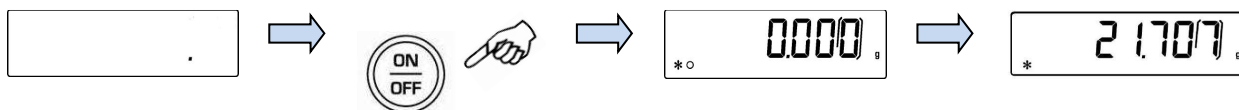

Press again **ON/OFF** button to return to “**STAND BY**” mode.


## 6. TARING

Place the container on the pan and press **O/T** button. “**O-t**” is displayed. As soon as the stability is reached the value of zero “**0.000**” will be displayed (weight of container is internally stored). Place the goods to be weighed in the container, the net weight will now be displayed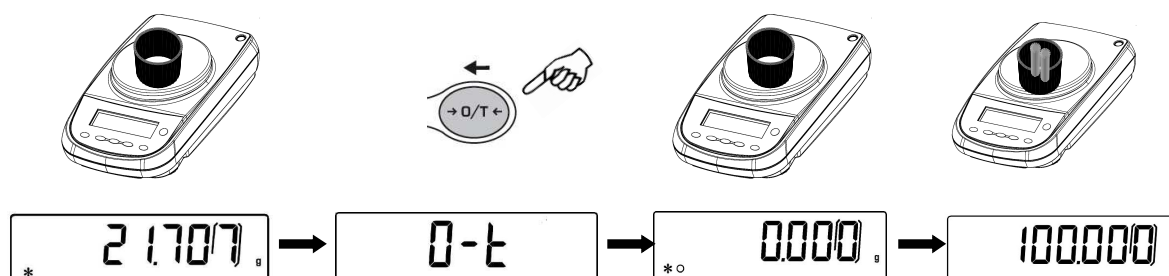


## 7. EXTERNAL CALIBRATION


Press the **CAL** button when no object is on the weighing pan; “**CAL**” will be displayed.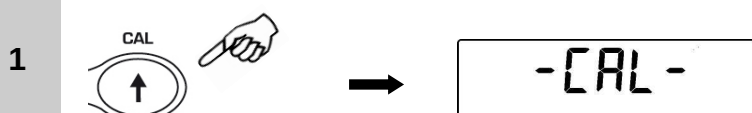

Wait until the value of required calibration weigh appears flashing, then put the required calibration weight carefully in the center of the weighing pan.

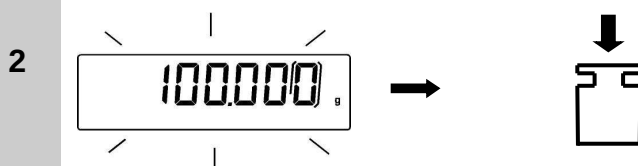
The display will stop flashing. After calibration is performed, the value of calibrated weight is displayed. Remove the calibration weight. Balance automatically returns to weighing mode and it is ready for weighing operations.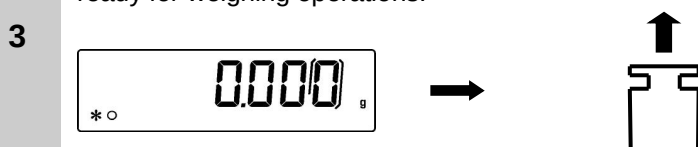

**NOTE:** if there is any interference during calibration process, an error message will be visualized.

## 8. OVERVIEW OF BALANCE PARAMETERS MENU

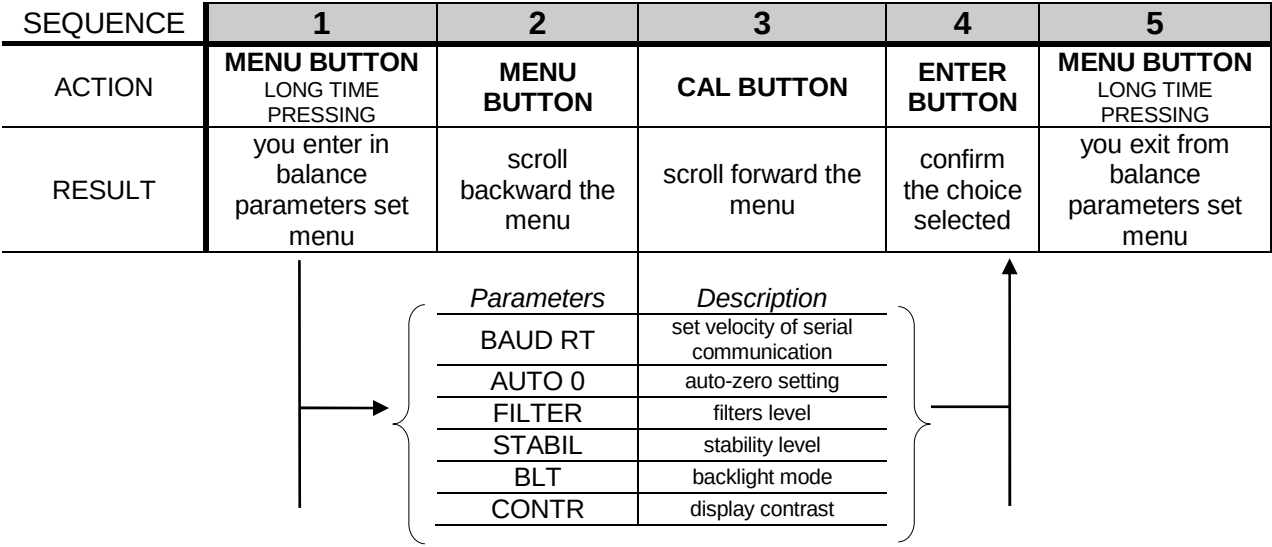
For a complete description of functions, it is warmly suggested to read the user manual of the balance

## 9. Quick guide to the use of the balance's programs

- To enter the menu of balance programs press the **MENU** button.
- Then use the **MENU** button to go to the next parameter, use the **CAL** button to return to previous parameter, use the **PRINT** button to confirm your selection.
- To escape from the menu press and keep pressed the **MENU** button until the acoustic sound is over.

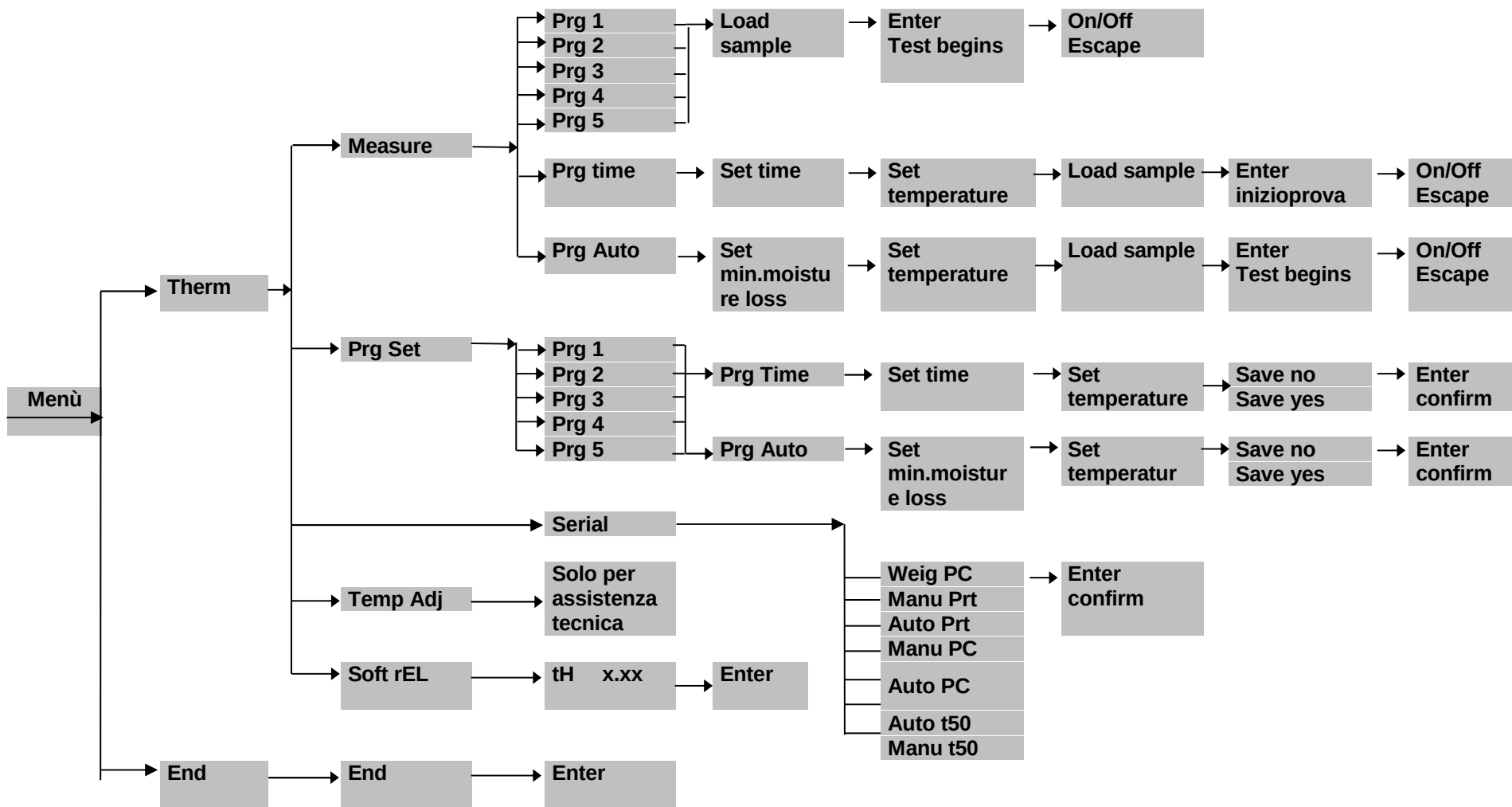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